


IndustrialTM Plankton

PBR 100L Installation Manual

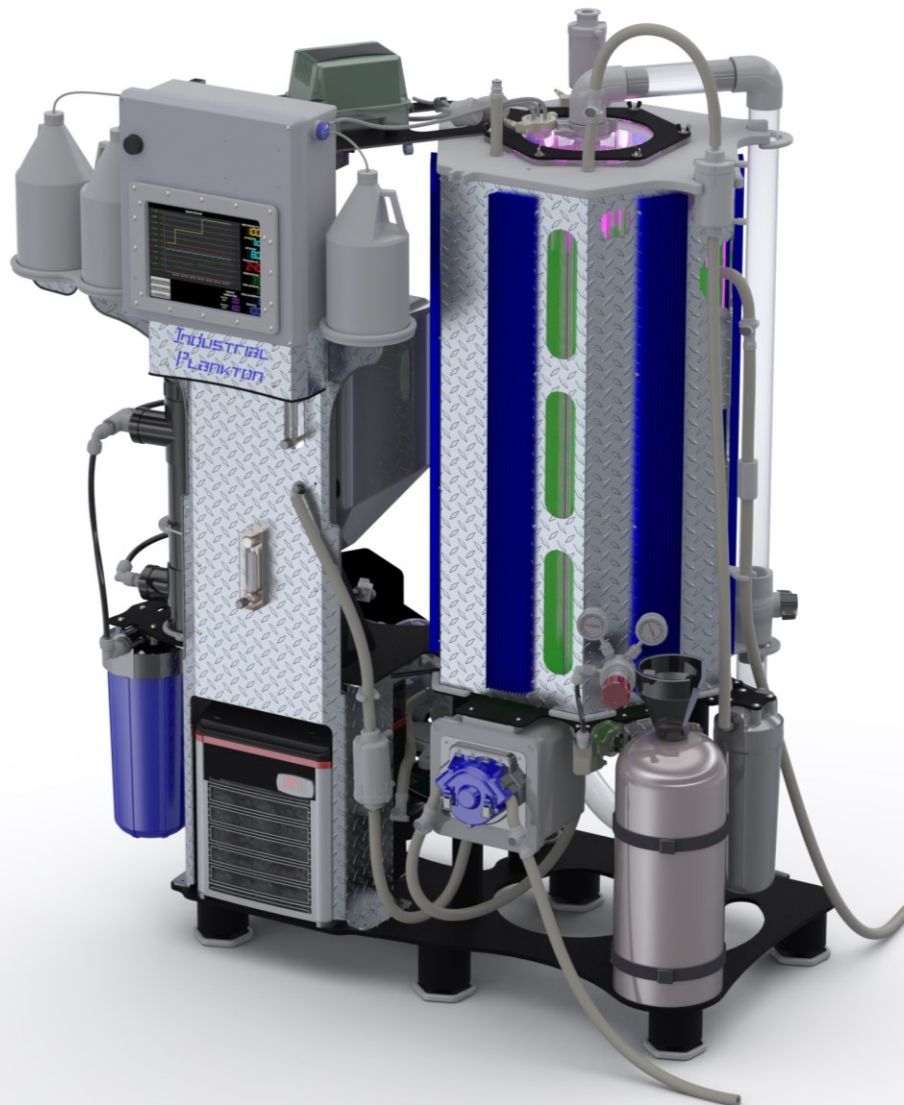

Warnings









1. **Never** use **Windex, acetone, ammonia,** or other solvent based cleaners on or near the PBR. The chemicals will weaken the acrylic and can cause catastrophic failure.
2. **Never** run Spray Pump without water in the PBR.
3. **Never** connect LED panels when the lights are powered. This damages the LEDs.
4. **Never** run Harvest Pump in reverse after inoculation.
5. **Never mix** cleaning fluids, it can result in serious injury or death.
6. **Never** drain the PBR without allowing air to replace draining fluid.
7. **Never** run the Chiller without water motion.
8. **Do not** drop the PBR when transporting it with a pallet jack or forklift.
9. **Do not** use sponges or paper towel to clean the tank. Only use soft cotton cloths to avoid scratching the acrylic.
10. **Do not** install PBR on soft ground or with uneven support.
11. **Do not** use the top port to add dangerous chemical without appropriate personal protection.









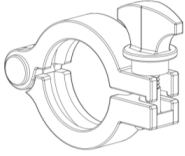
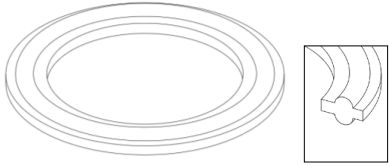
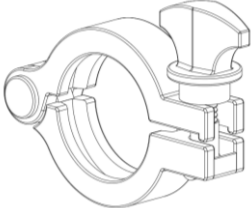
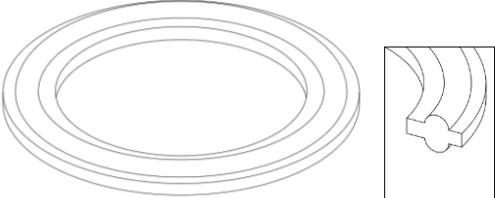


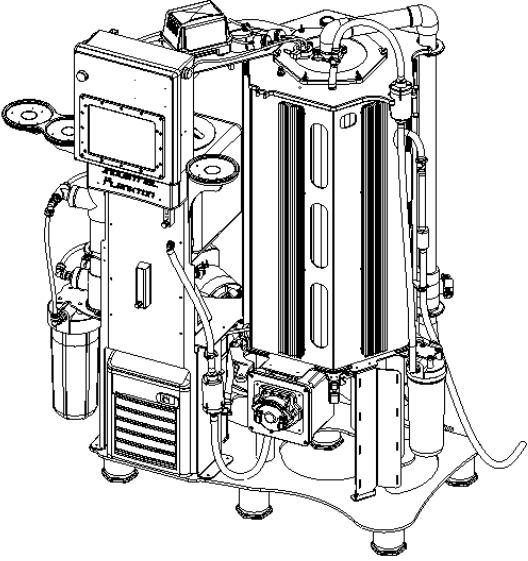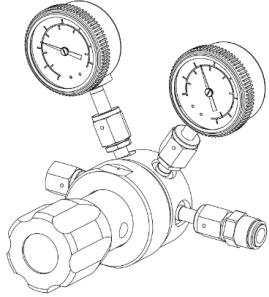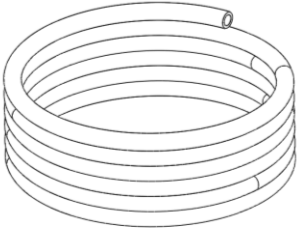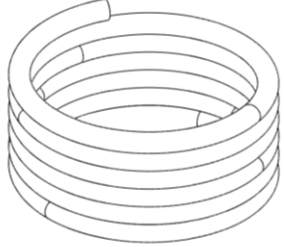Table of Contents

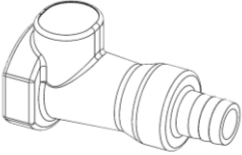PBR 100L Installation Manual.....	1
Warnings.....	2
Table of Contents	3
Contents of Crate	4
Spare Parts Kit	6
Abbreviations Used	6
Uncrating	7
PBR 100L Components (Front)	8
PBR 100L Components (Back).....	9
Place the PBR.....	10
Water Flow	11
Water Connections.....	11
Drainage	15
Harvest Lines	15
Temperature Control.....	17
CO ₂	18
Top Reassembly.....	19
Power.....	19
Wi-Fi	20
Advanced Wi-Fi Setup	20

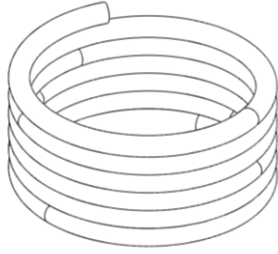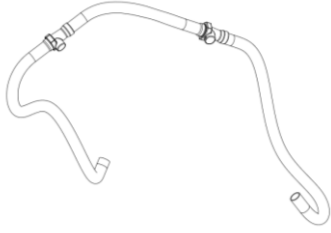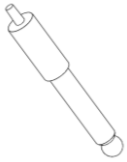Contents of Crate

 <p>PBR 100L R1006 x 5ft</p>	<p>P1099 x 1</p>  <p>CO₂ Regulator (with 3/8" JG fitting)</p>
 <p>CO₂ Tubing: 3/8" Black Poly Tubing B1191 x 1</p>	<p>R1082 x 7.6m (25ft)</p>  <p>Water Inlet Tubing: 1/2" Black Poly Tubing</p>
 <p>Hose Clamp Pliers (for constant tension hose clamps)</p>	<p>B1235 x 1</p>  <p>1/2" mNPT x 1/2" JG Fitting (attached to R1082)</p>
<p>B1195 x 2</p>  <p>Female 3/8" Quick Disconnects (3/8" fQDC) (For inoculant carboys)</p>	<p>B1034 x 3</p>  <p>Stir Bars</p>

<p>R1001 x 25ft</p>  <p>Drain tubing: 3/4" ID vinyl hose</p>	<p>S1052, S1053</p>  <p>Automated Harvest Line with Extension</p>
<p>B1289 x 1</p>  <p>pH Probe (Ensure fluid is in cap)</p>	<p>B1429 x 1</p>  <p>140 mL Syringe</p>
<p>P1037 x 1</p>  <p>Pleated 0.2 µm Filter</p>	 <p>10" 1 µm Filter</p>
<p>P1042 x 1</p>  <p>Air Vent Filter</p>	 <p>UV Sterilizer Bulb</p>
<p>B1163 x 1</p>  <p>1.5" Triclamp</p>	<p>B1037 x 1</p>  <p>1.5" EPDM Triclamp Gasket</p>
<p>B1164 x 1</p>  <p>2" Triclamp</p>	<p>B1206 x 1</p>  <p>2" EPDM Triclamp Gasket</p>

Spare Parts Kit

PN	Description	QTY
B1364	Silicone Lubricant	10 mL
B1037	EPDM Sanitary Flange Gasket 1.5"	1
B1206	EPDM Sanitary Flange Gasket 2"	1
B1094	1/4"-20 wingnuts for lid	4
B1089	Neoprene roofing washers 5/16" ID	4
B1103	#10-32, 3/4" button head screw	2
B1030	Constant tension tube clamp for 1/2" ID crush resistant silicone	3
B1217	Spring clamps for harvest pump hose (3/8" ID)	3
B1311	Constant tension spring clamp for nutrient tubing (3/16" ID)	3
B1011	Locking clip for 3/8" OD push-to-connect tubing	3
B1313	Locking clip for 1/2" OD push-to-connect line	3

Abbreviations Used

Abbreviation	Description
PBR	PhotoBioReactor
PLC	Programming Logic Controller
QDC	Autoclavable Quick DisConnect fitting
mQDC	Male autoclavable Quick DisConnect fitting
FQDC	Female autoclavable Quick DisConnect fitting
NPT	National Pipe Thread
F/2	Guillard's F/2 algae growth media
CIP	Clean-in-place
GFI	Ground Fault Interrupter
thio	Sodium thiosulphate

Uncrating

Required Tools

- PBR½” Socket wrench or crescent wrench
- Pallet jack or equivalent
- Compressed gas CO₂ cylinder
- Drill or changeable bit screwdriver


1. Unscrew Drawbridge Wall of the crate.




2. Remove loose packages from beside PBR and locate #2 Robertson (square) driver.




3. Unscrew hardware securing the PBR feet to the crate floor

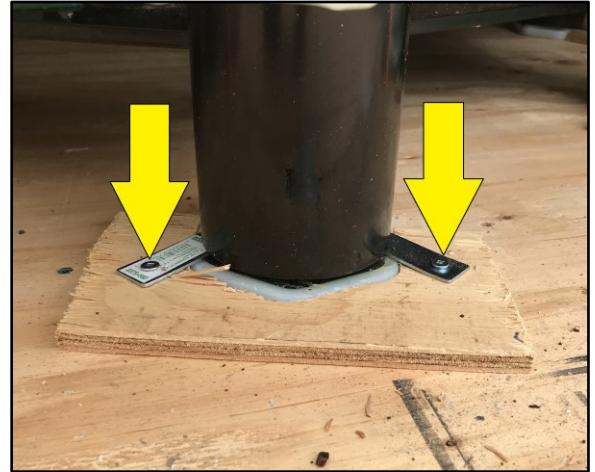


4. Use a pallet jack or forklift to remove the PBR from the crate. **Ensure not** to scratch the powder coated base, to avoid corrosion.

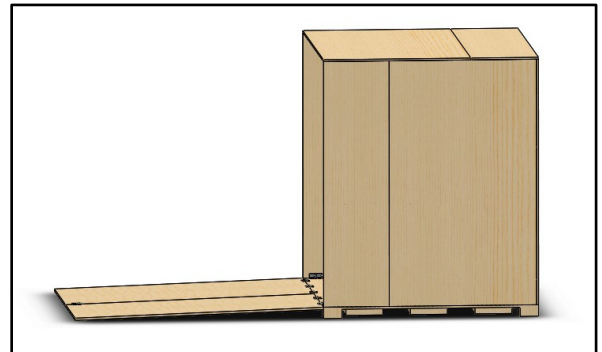

PBR 100L Components (Front)

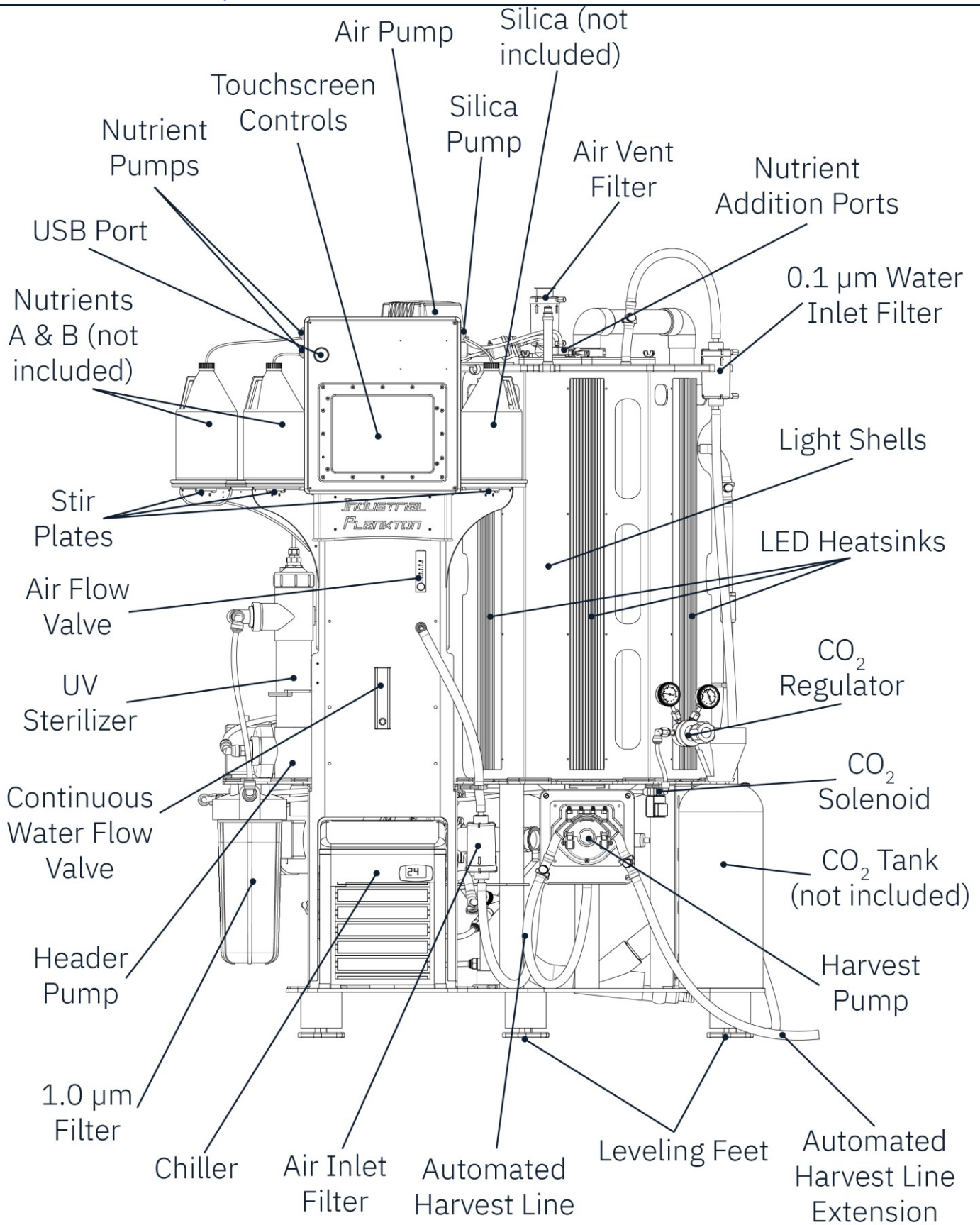

PBR 100L Components (Back)

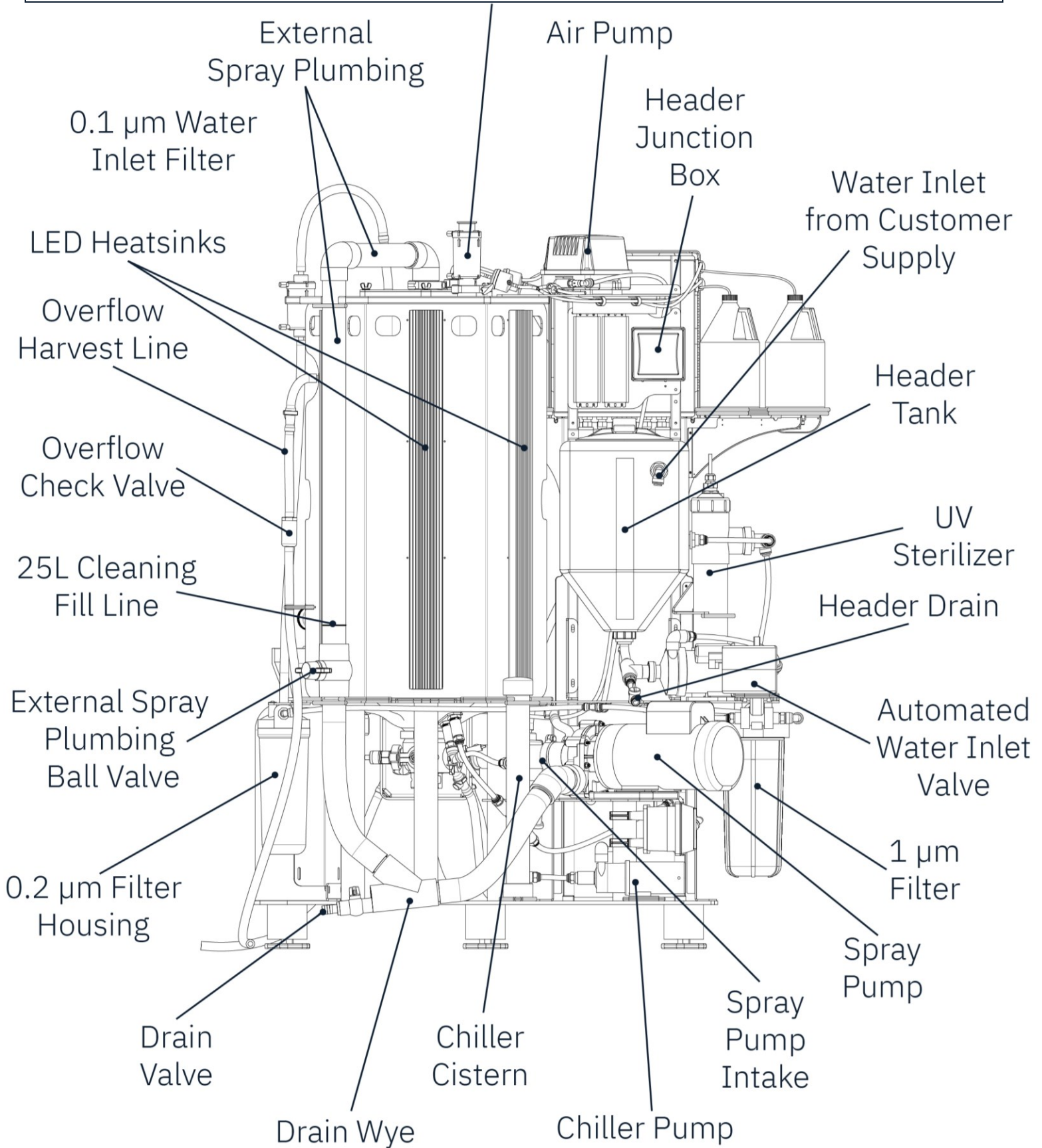


Place the PBR

Leveling and evenly supporting the PBR is critical for spray system functionality and lifetime of the tank.


Required tools:

- Pallet jack or similar
- Spirit/bubble level
- Paper
- IP Foot Wrench

1. Use spirit on the lower steel plate to measure level. Ensure to check level in several different orientations.

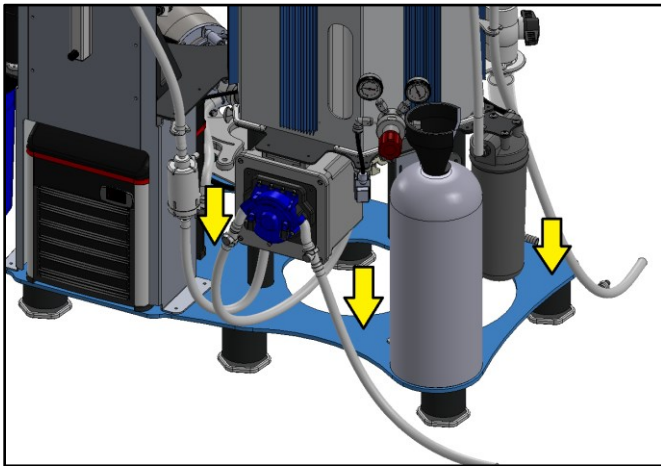


2. Rotate the leveling feet to adjust to the PBR to level.

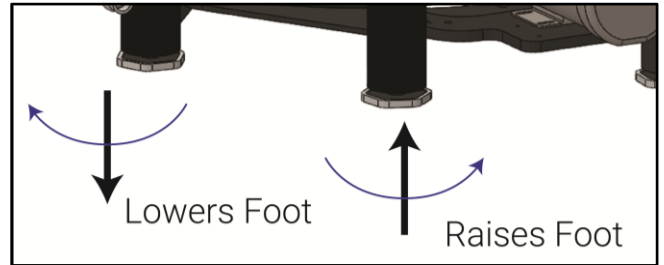

3. Once level, ensure each foot is supporting the PBR evenly. Slide a thin sheet of paper under each foot to ensure it is tight to the ground – paper should not slide between the foot and the floor.



Water Flow

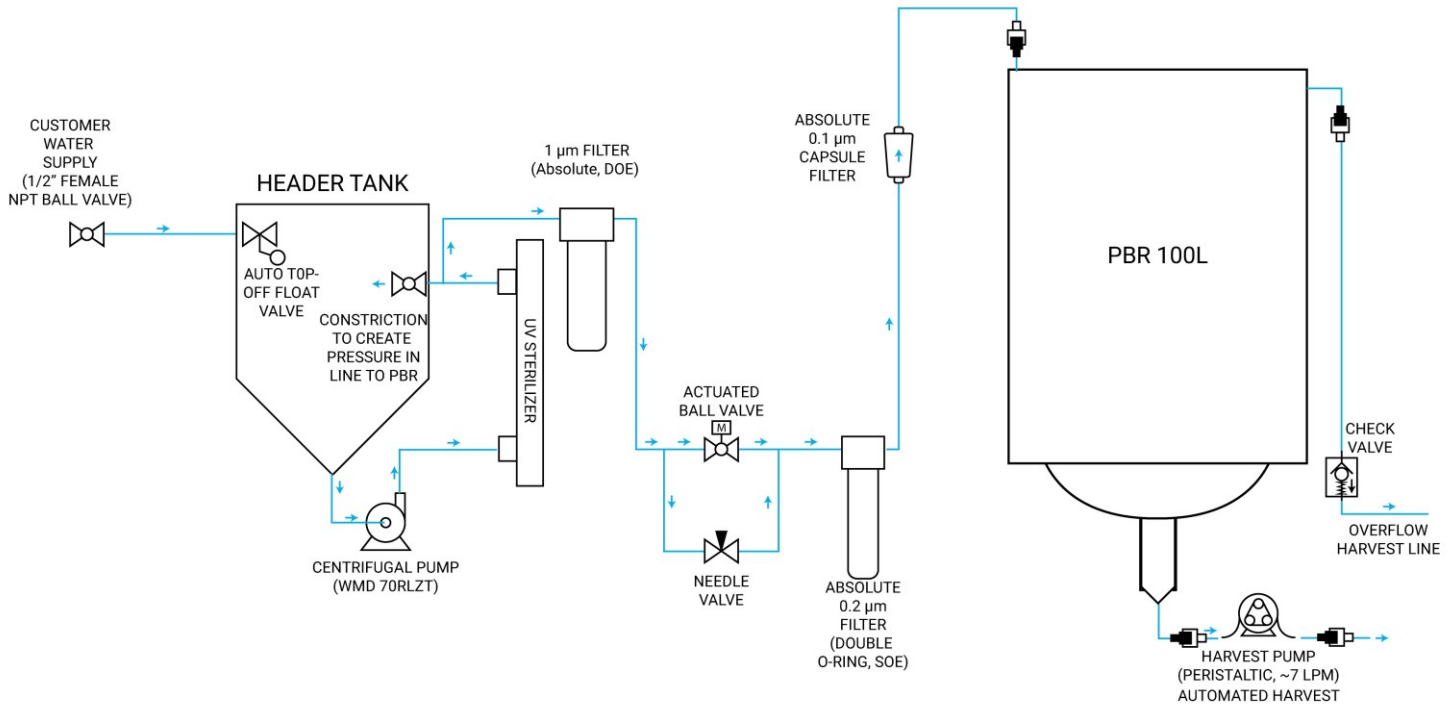

Water Connections


Ensure that push-to-connect tubing is fully inserted into fittings to create a proper seal. If there is a leak, push tube in further. To remove a tube, depress the dark grey ring toward fitting body and pull on tube.

1. Tighten supplied 1/2" mNPT by 1/2" OD push to connect adapter (B1235) into the facility's water supply connection. Use plumbers (PTFE) tape to seal threads.
2. Ensure end of supplied 1/2" OD black poly tubing (R1082) is round and cut square, then insert into B1235.

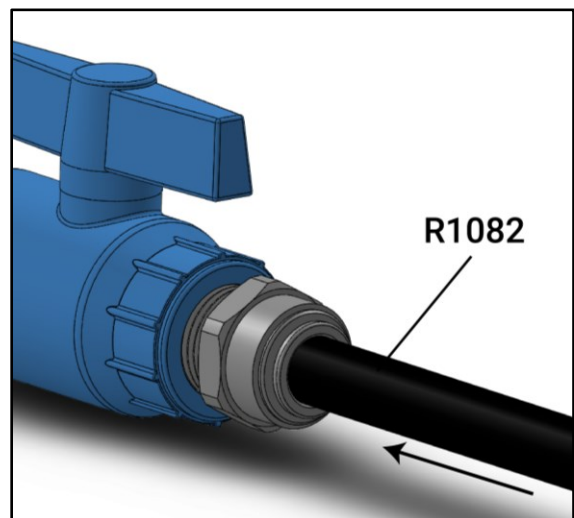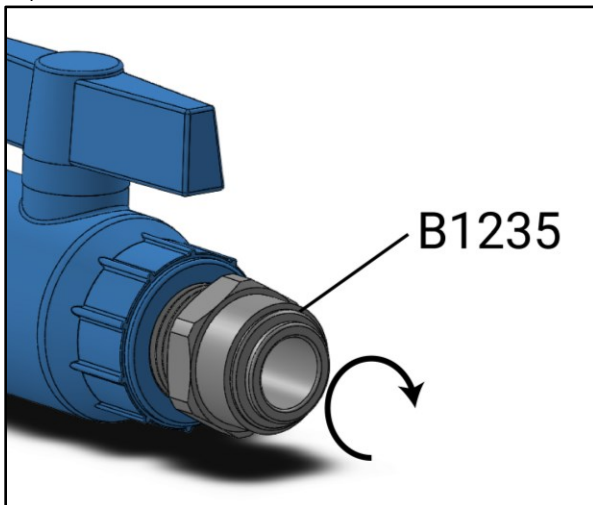


Industrial[™] Plankton


3. Unroll and route R1082 connected to water supply (from step 1) to the float valve bulkhead on the Header Tank.

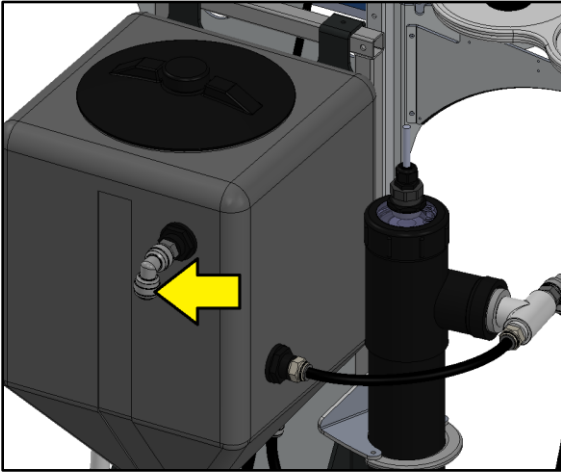


4. Loosen UV cord grip on Header Tank.

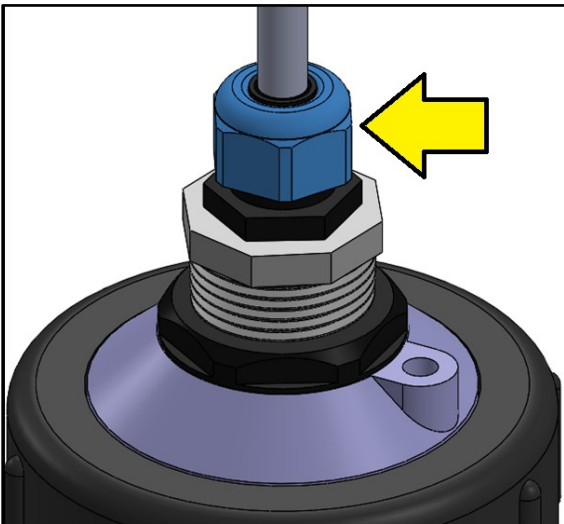


5. Unscrew UV cord fitting and remove plug from UV sterilizer.

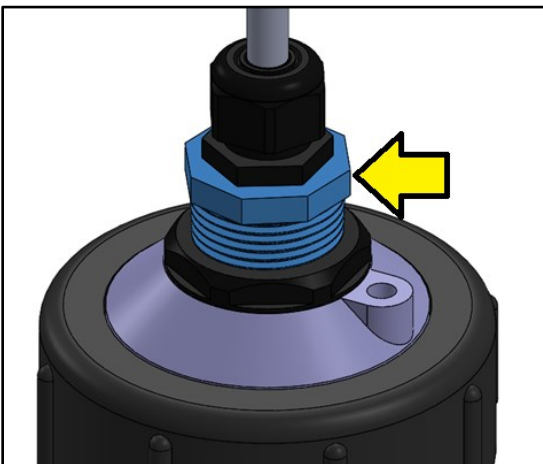


6. Insert UV bulb into Quartz sleeve and connect the 4-pin connector to the bulb. Do not touch bulb's glass with bare hands or drop the bulb.

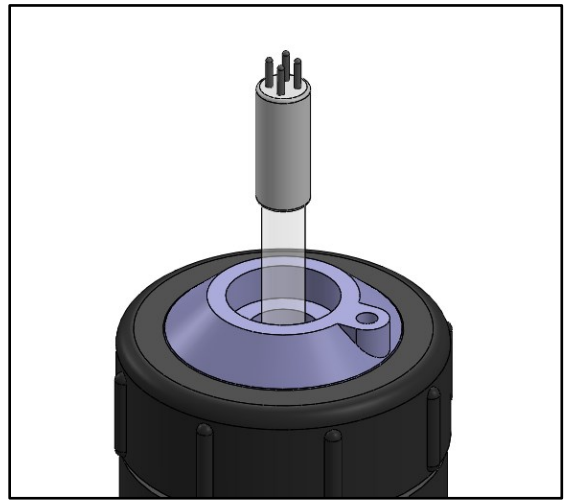


7. Tighten UV cord fitting into UV Sterilizer. If water enters the quartz sleeve it will damage the bulb and ballast.

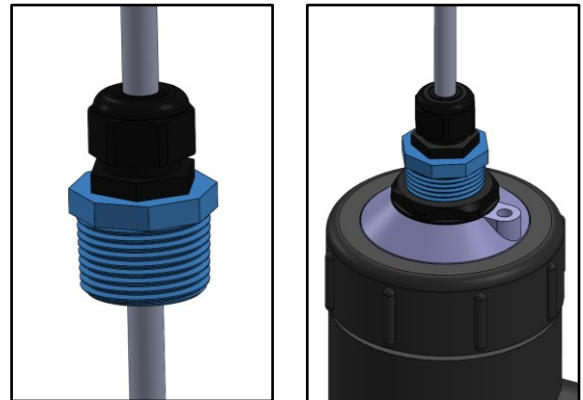


8. Tighten cord grip.

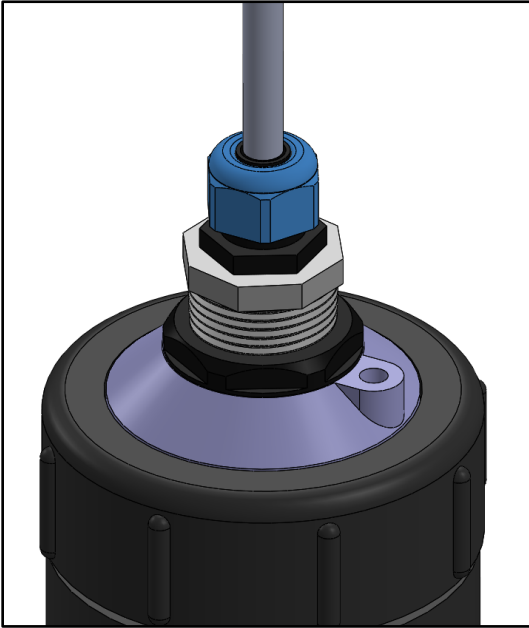

9. Loosen canister on 1 µm water filter housing.

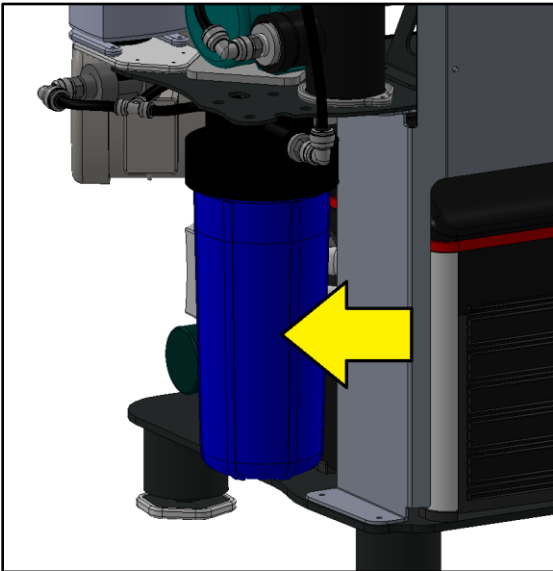


10. Place 1 µm water filter into bottom of filter housing ensuring it is centred.

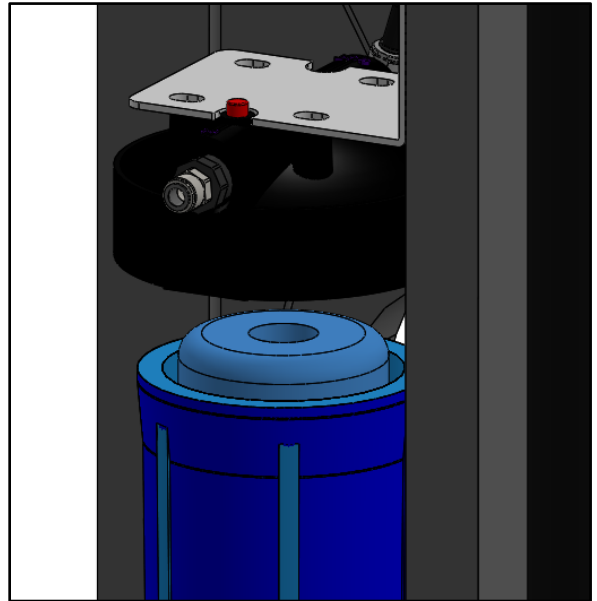


11. Keep filter centred and screw canister back into top.

12. Unscrew 0.2 µm filter housing canister on PBR and check O-ring.

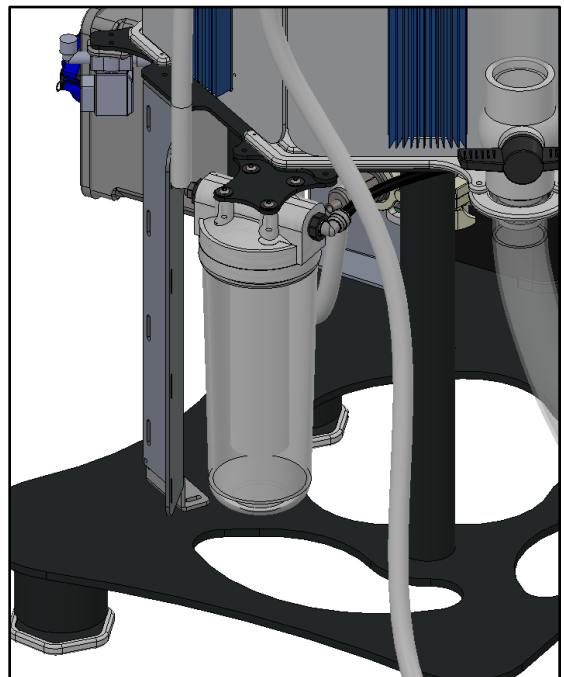


Industrial™ Plankton

13. Place filter membrane into canister below filter top (double O-ring up)

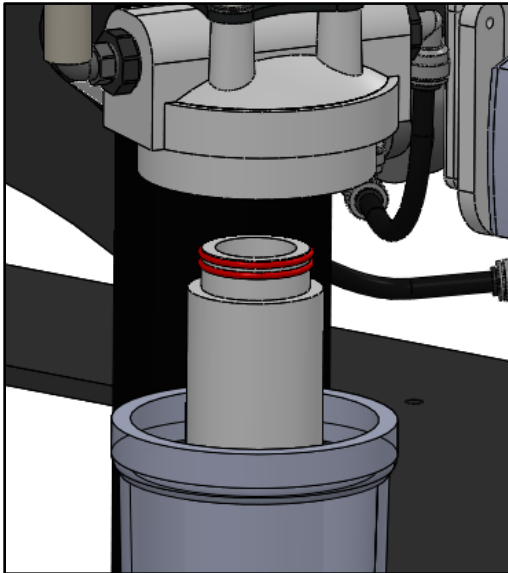

14. Fully insert filter membrane into top

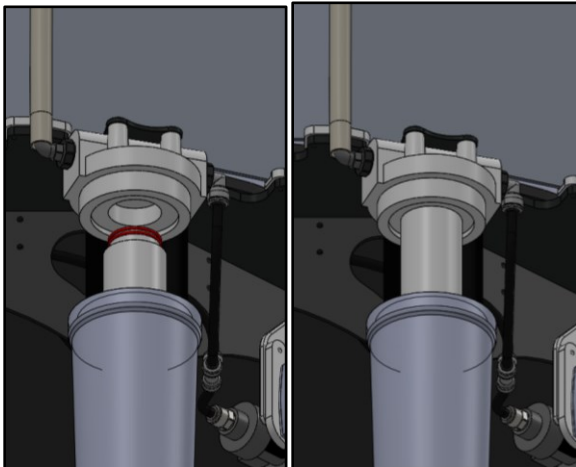


15. Ensure O-ring is seated in canister to avoid shearing, then tighten closed.

16. Turn on water supply

17. Check for leaks in all water lines


18. Bleed air from 1 μm (red button) and 0.1 μm filters (top side-barb).

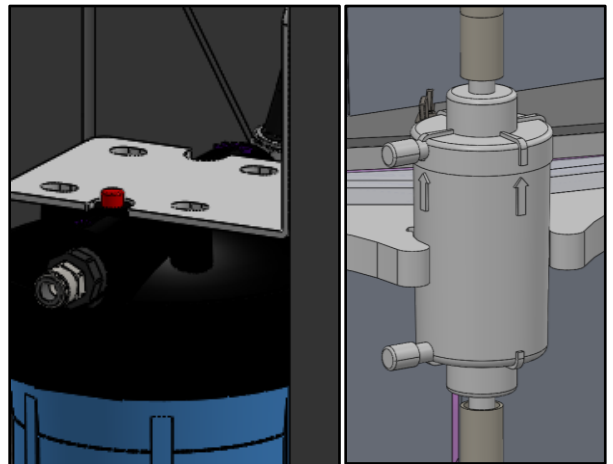

Drainage

The PBR is designed to gravity drain. If a floor drain is unavailable, the harvest pump can be used to drain the PBR. Refer to tank “Harvest: Manual” description in the user manual for details.

1. Attach **R1001** (¾” ID vinyl hose) to drain valve
2. Route **R1001** to the floor drain

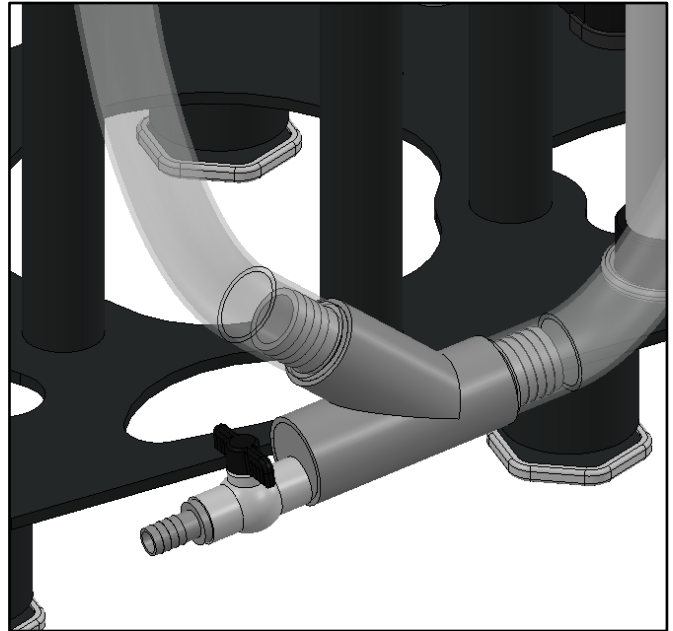

Harvest Lines


Never submerge Overflow Harvest Line or run uphill. It relieves excess air pressure inside the tank to stabilize volume readings and prevent top seal from leaking during spray cleaning.

Never connect the Overflow Harvest Line to waterways with livestock when cleaning. Cleaning agents must go to waste.

The Overflow Check Valve prevents contaminated air from entering the PBR when harvesting/draining. **Ensure** air and water can flow freely out of the PBR through it.


1. Route the Overflow Harvest Line to algae collection tank (or where algae is desired) when culturing in continuous mode. Or, leave directed to waste for spray cleaning. Any tubing or vessel attached to this line must be

able to vent to ambient pressure.

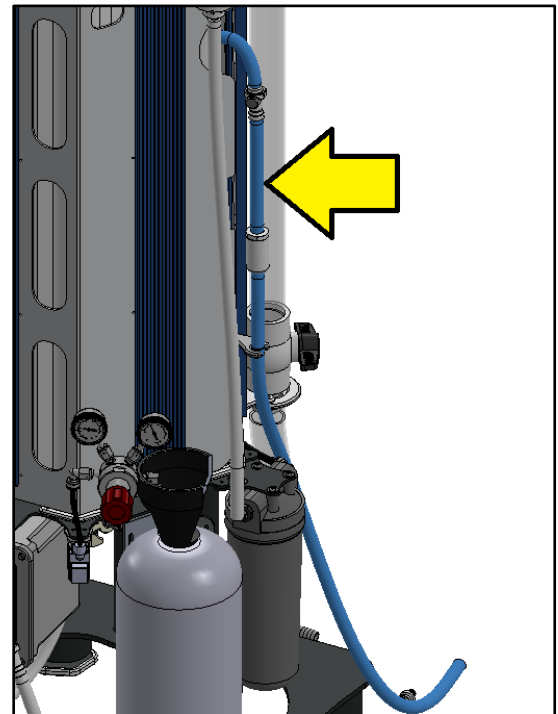


INDUSTRIAL™ PLANKTON


2. Attach Automated Harvest Line to bottom barb on Manifold and adjust tube clamp to secure it. Ensure mQDCs are closer to the PBR than fQDCs.

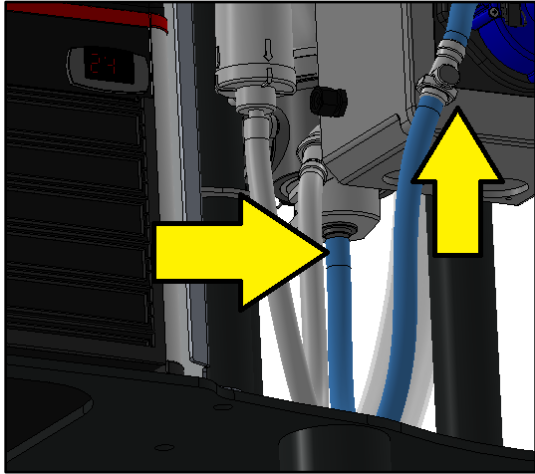


3. Unclamp and open Harvest Pump top cover.

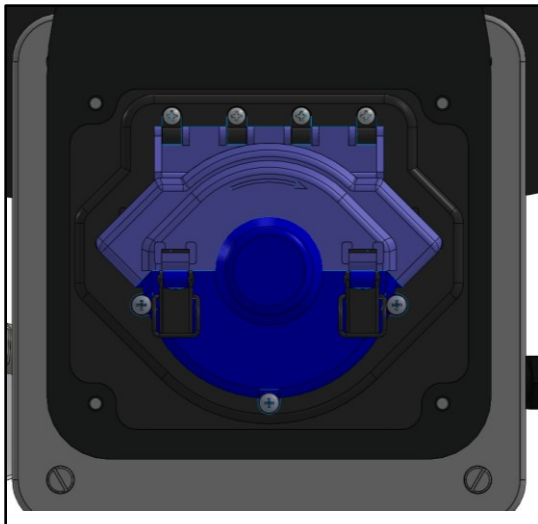


4. Apply **B1364** (silicone lubricant) to the Automated Harvest Line section between QDCs.

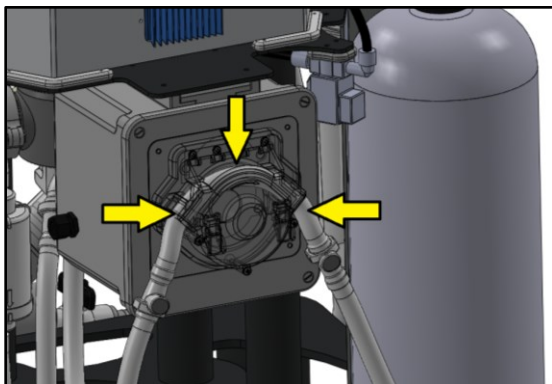

5. Place this section through the Harvest Pump's rollers, ensuring to press the tubing far back into the rollers. Ensure that the tube is pressed back, past the protruding tabs, to avoid damage to the tube and pump.

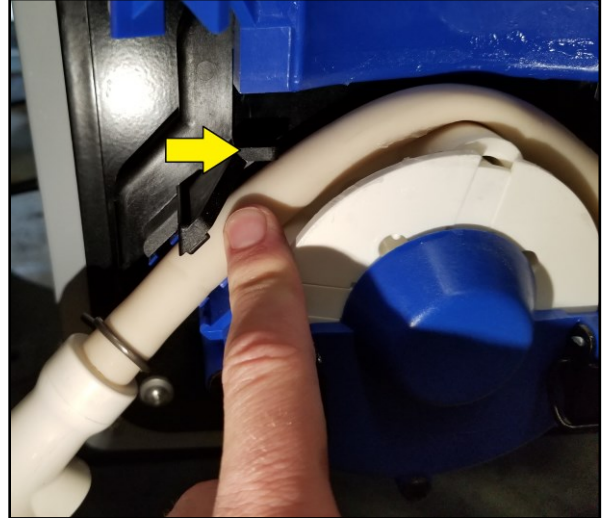


6. Press top cover down and clamp shut.

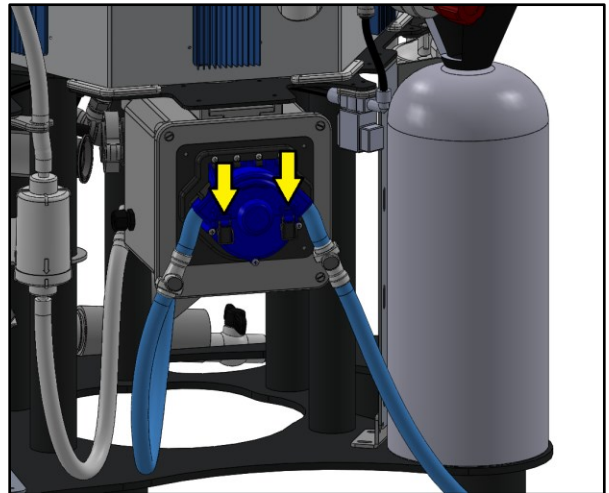


Temperature Control


- Warnings:
 - Never run chiller without water motion
 - Carry chiller by handles
 - Top up chiller cistern monthly or as required

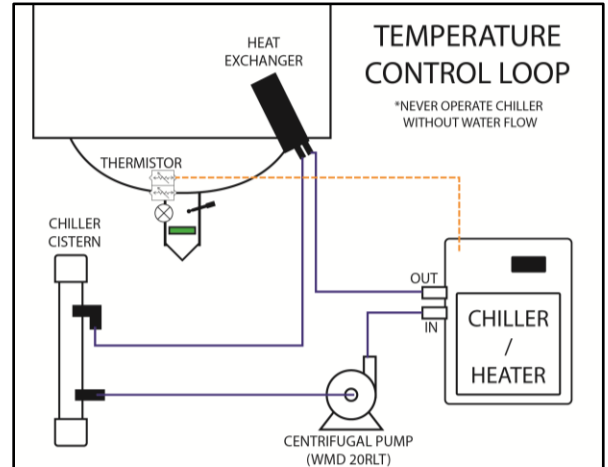

1. Remove Chiller Cistern Cap

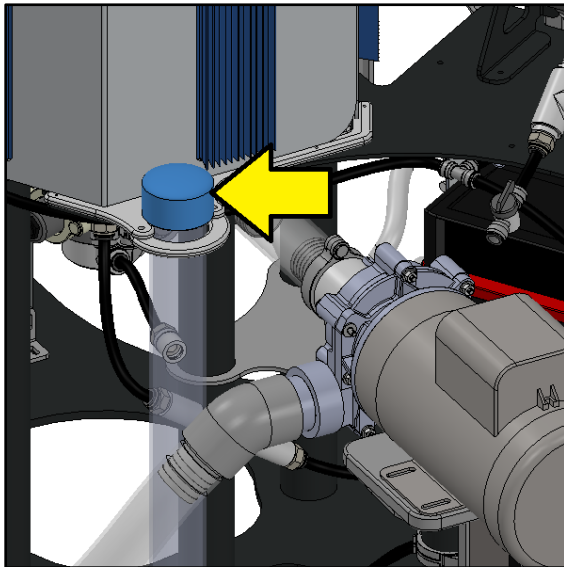


2. Fill cistern to Fill Line

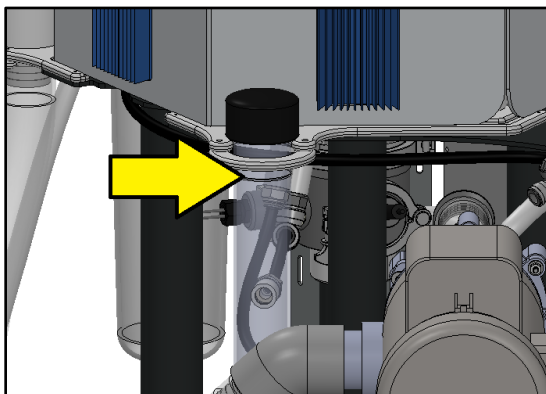


3. Energize the Chiller Unit (the junction box powers both the chiller and chiller pump) and top up cistern (air will be trapped in the lines). It may be necessary repeat several

times.

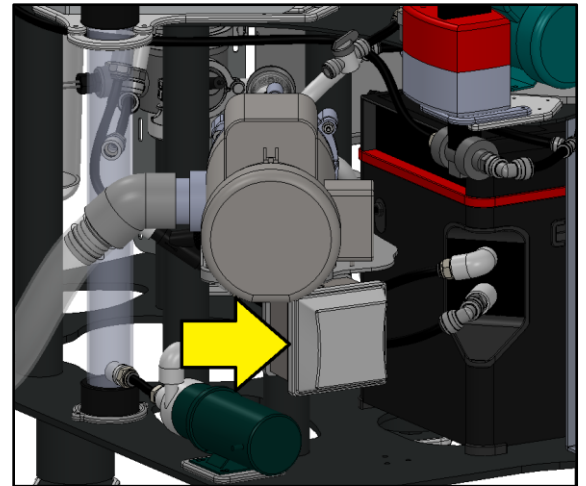


4. Replace cistern cap to avoid evaporation and check for leaks

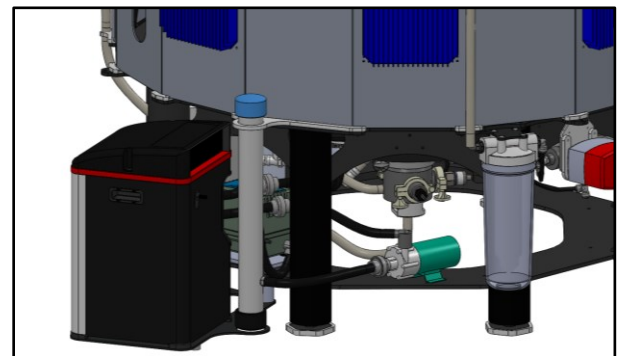


5. Turn on chiller. For instructions refer to [Teco TK 500 / 1000 / 2000 manual](#)

CO₂


1. Place CO₂ cylinder (**not included**) in holder and tighten straps to secure in place.

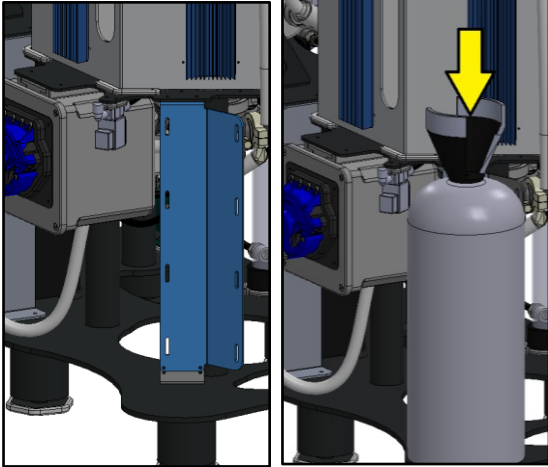

2. Attach **P1099 (regulator)** to CO₂ cylinder (check gasket for debris).

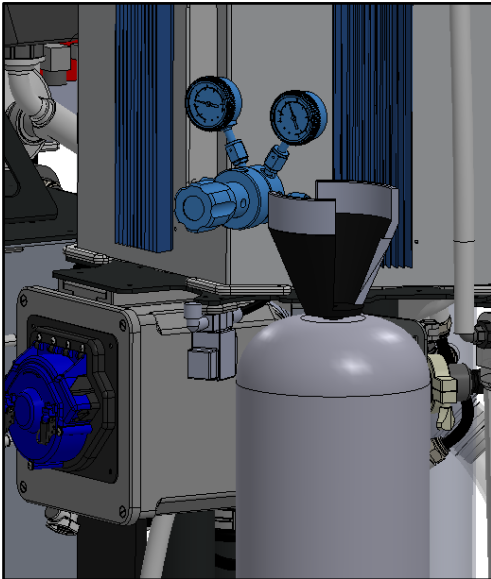


3. Connect regulator to CO₂ valve with **R1006 (3/8" black tubing)**

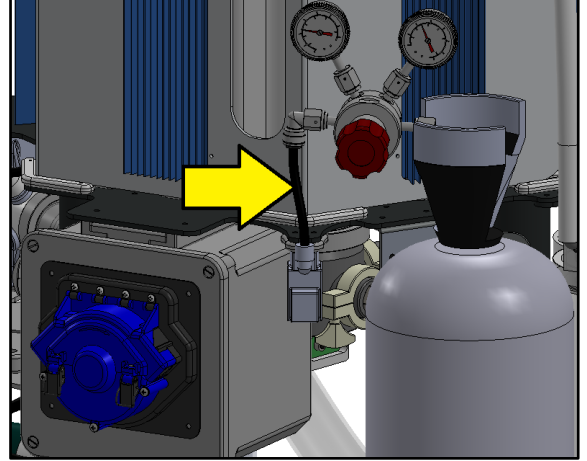


4. Open CO₂ cylinder valve fully.
5. Check regulator is set to 15 psi (pull out red knob to adjust and push back in to set).

Top Reassembly


1. Attach Air Vent Filter (P1042, KA3PFRP1) to triclamp with gasket (B1037) and clamp (B1163) on lid with flow direction going out of the tank.

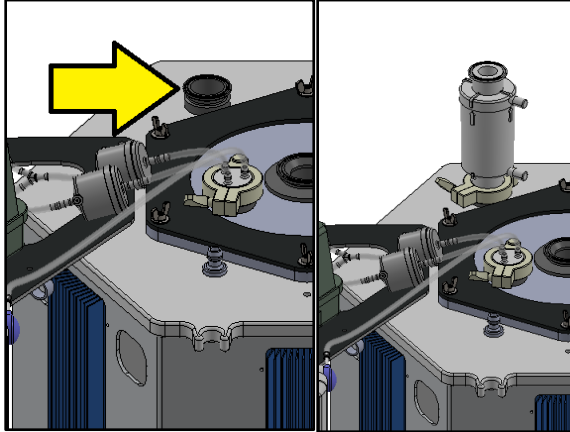

2. Connect external spray post using 2" triclamp (B1164) and gasket (B1206)

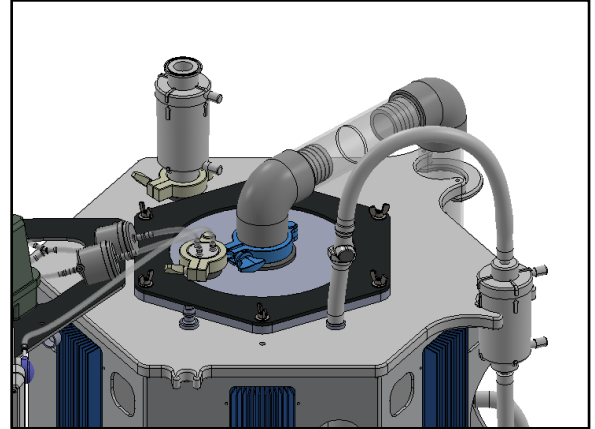

3. Connect top air quick disconnect air tube

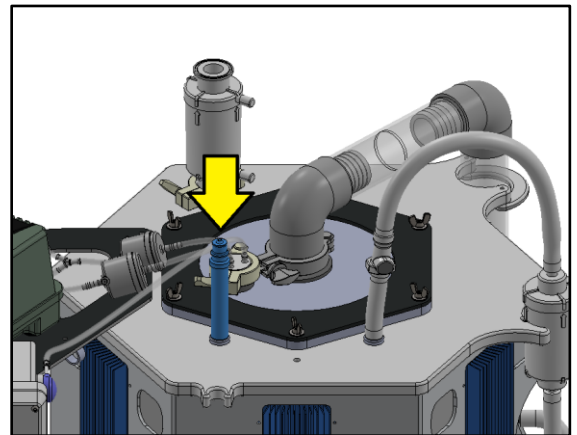


Power

Plug in the reactor and it will automatically boot into the algae program.

Wi-Fi

To connect to the PBR remotely, Wi-Fi internet must be available, and your remote access subscription must be current.

1. Go to Advanced Settings.

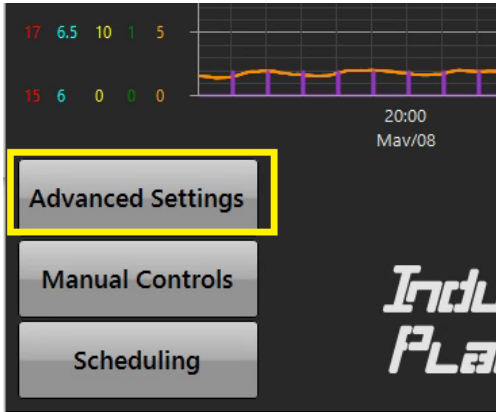

2. Press Continue.




3. From Advanced Settings select Wi-Fi network from list and enter the password.




4. If step 3 doesn't connect, continue to [Advanced Wi-Fi Setup](#).

Advanced Wi-Fi Setup

1. Go to Advanced Settings

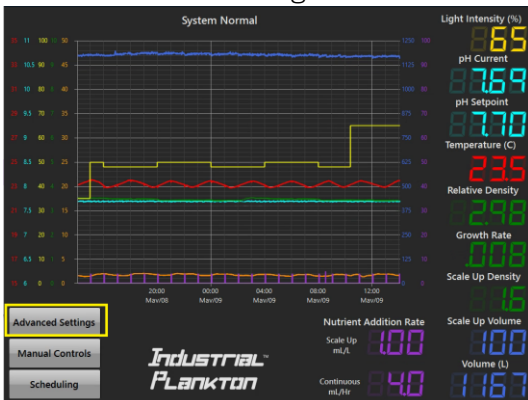


2. Press Continue.

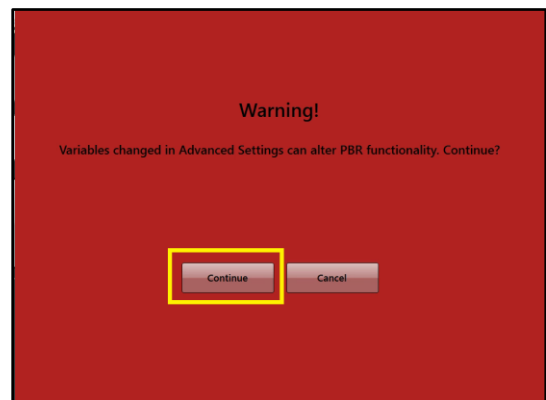

3. Press Exit Program.




4. Open the on-screen keyboard (in lower right menu on windows login screen).

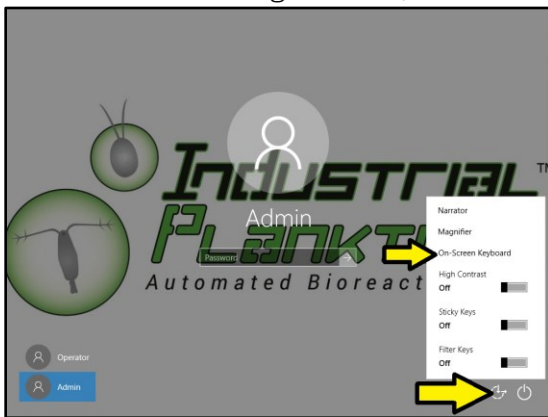


5. Log into Admin (password: dansarki).

6. Double click the Unlock computer icon on the desktop (this will restart the computer).

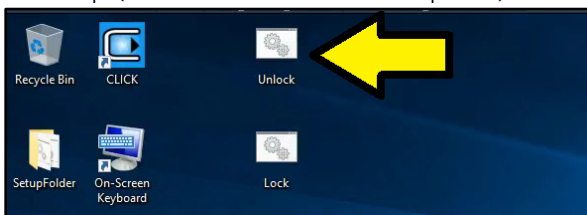


7. Repeat steps 1-5.


8. Open On-Screen Keyboard (only required if network is password protected).

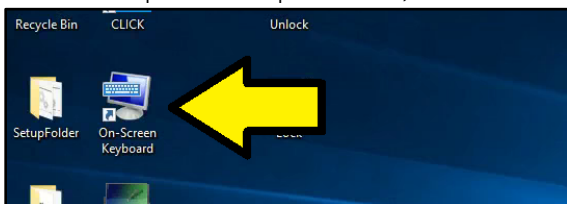

9. Drag the keyboard to the left of the screen and select Wi-Fi Logo in right system tray.

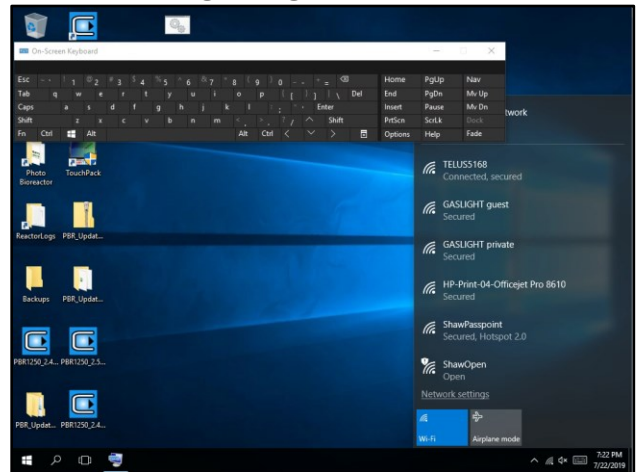

10. Select network and enter password.

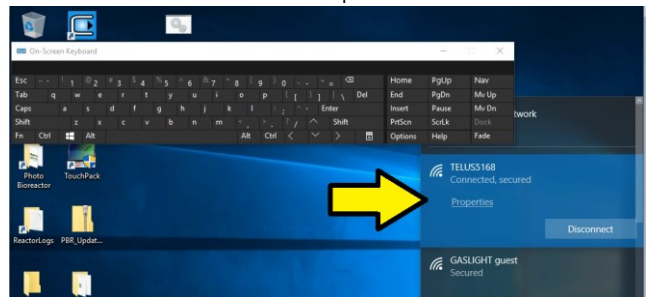

11. Double click "lock computer" icon on desktop.



12. The PBR will restart and launch the control system upon reboot.